



1. Turning off the system.


- a. Exit Zen Black – a pop-up window will appear showing which lasers are on. Click all lasers that are on to the off position.
- b. Shut down Acquisition Computer.
- c. Wait about 15 min for the Fan on the HeNe Laser to stop running, then turn off the Main Power Switches in the reverse order to that used during startup (Components, System PC, then Main Switch (see image below)).

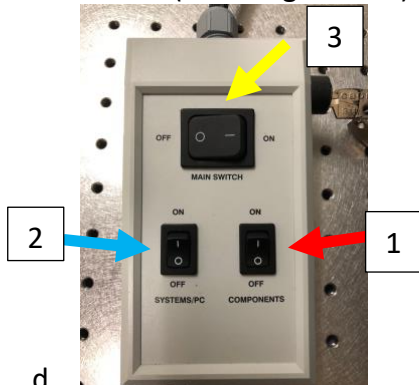

- d.
- e. Retrieve data from Analysis Computer
 - i. Image files are located in the DATA (D:) drive,

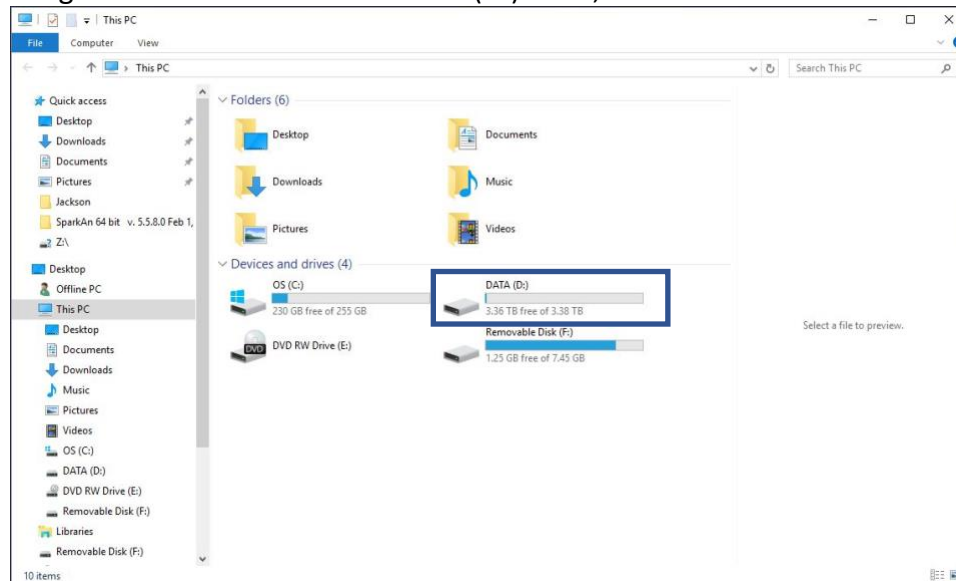

- ii.
- iii. In the path: This PC > DATA (D:) > SWAP > Zeiss > Your Folder as shown below

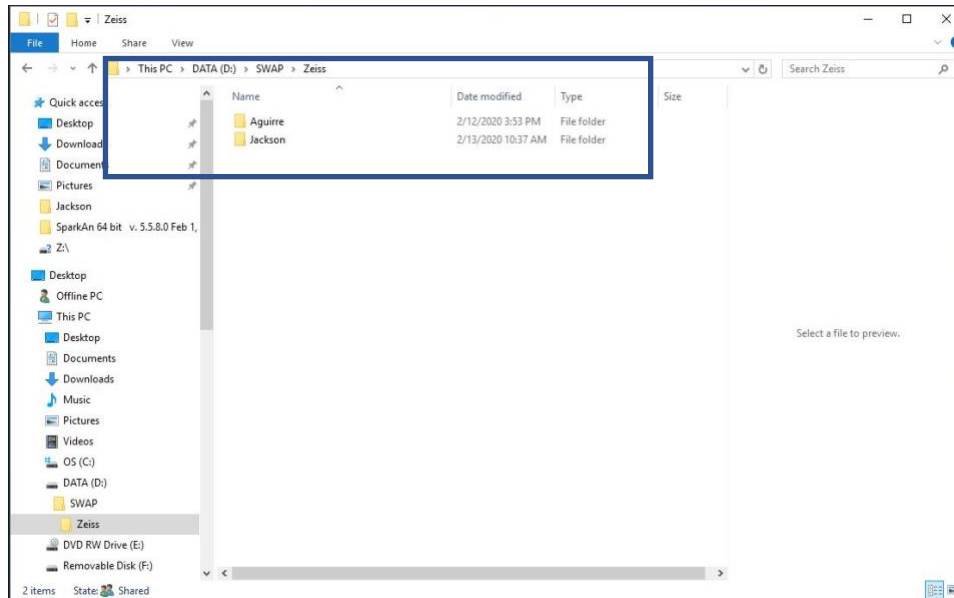

- iv.
- v. Where Your Folder is the name of the folder for your lab (see Dr. Jackson if you need help).
- vi. Either use a large capacity Flash Drive (64 GB or larger), external USB drive or
- vii. Send files to your One Drive, Google Drive, or other cloud storage (this is very slow).
- f. When files have transferred, shut down Analysis Computer.